



13th International mRNA Health Conference @Berlin

Novel Ionizable Lipids **for in vivo liver, ex vivo CAR knock-in and in vivo liver-detargeting**

Shinichi HASHIMOTO

Research Manager, Head of LNP Formulation Discovery

Bio Science & Engineering Lab.
FUJIFILM Corporation

11-13 Nov 2025



FUJIFILM
Value from Innovation

Forward Looking Statements and Regulatory Matters

This presentation contains certain statements which constitute “forward-looking statements”. These forward-looking statements may be identified by words such as ‘believes’, ‘expects’, ‘anticipates’, ‘projects’, ‘intends’, ‘should’, ‘seeks’, ‘estimates’, ‘future’ or similar expressions or by discussion of, among other things, strategy, goals, plans or intentions. The forward-looking statements involve risks and uncertainties that could cause actual business, financial, and technology, clinical and regulatory development results to differ materially from those expressed in the forward-looking statements. Many of these risks and uncertainties relate to factors that are beyond Fujifilm’s abilities to control or estimate precisely, such as future market conditions, the behaviors of other market participants, the technological success of Fujifilm’s preclinical- and clinical-stage programs, regulatory authorization or approval of Fujifilm’s product candidates, and other business effects, including the effects of industry, economic or political conditions, and therefore undue reliance should not be placed on such statements. Examples of forward-looking statements in this presentation include, but are not limited to, statements regarding the market for LNP-encapsulated drugs and biologics and the potential of Fujifilm’s LNP technology to result in one or more competitive products that are authorized or approved by applicable regulatory agencies in one or more countries. Actual results may differ materially from those in the forward-looking statements.

This presentation contains statements related to the biological, chemical, medical, and related characteristics of product candidates under development by Fujifilm and the commercial promotion, distribution, and sale of which will require authorization or approval from regulatory agencies on a country-by-country basis. Positive results of preclinical experiments in nonhuman laboratory models is no guarantee of such results in clinical studies in humans. None of the product candidates described in this presentation have been authorized or approved by any regulatory agencies; and nothing contained in this presentation should be regarded as the promotion or marketing of any such product candidates or any review or decision by any such regulatory agencies as to the safety or effectiveness of such product candidates, or whether such product candidates will be authorized or approved by any such regulatory agencies.

FUJIFILM Ionizable Lipids | Design concept and list

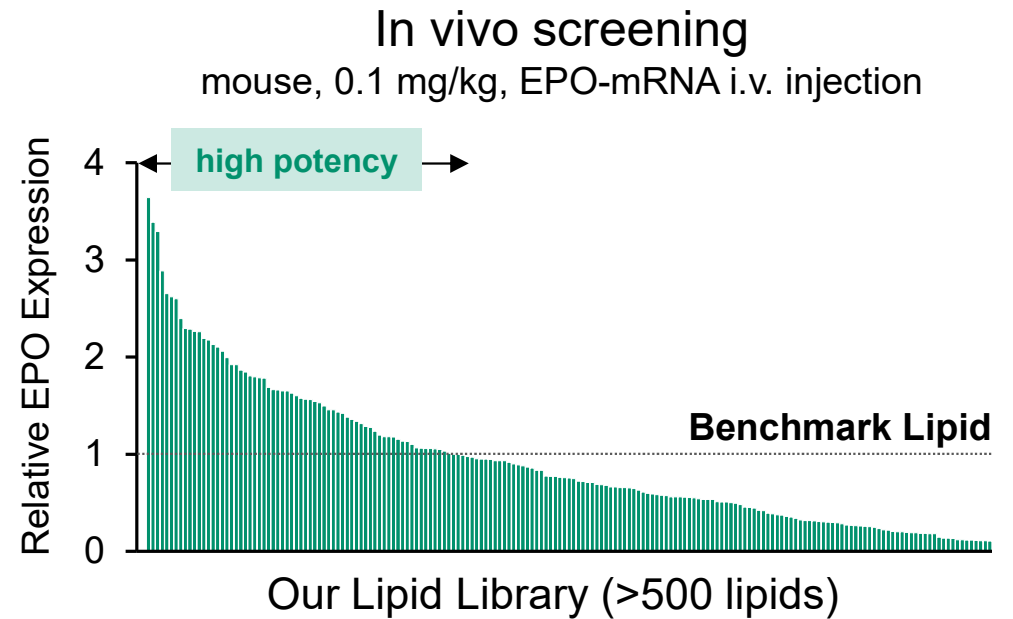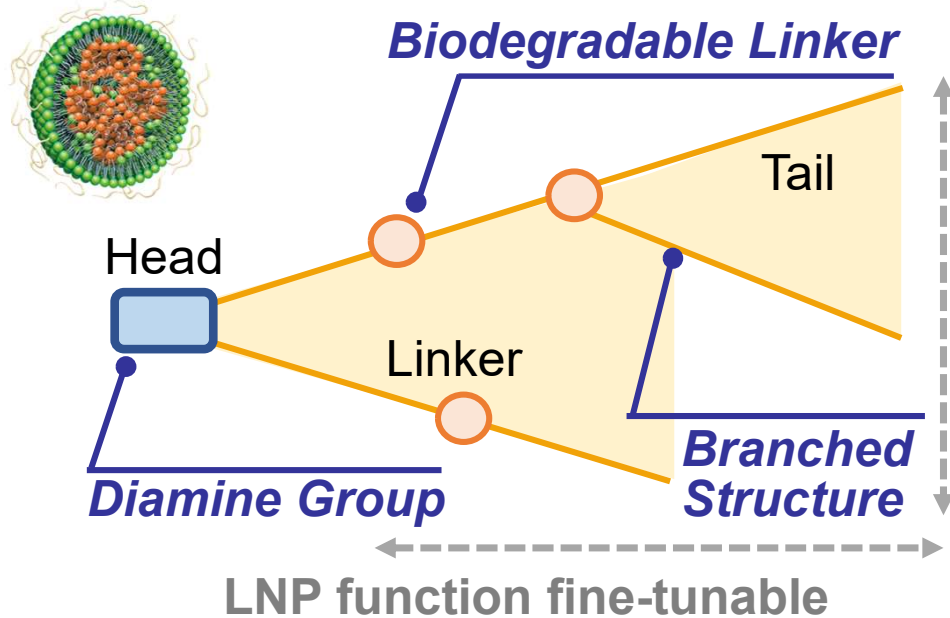

In vivo mRNA-LNP therapy

i.m.(vaccine)

FL-0445

GMP

i.v. liver

FL-1245T

New

extra-hepatic

FL-1779T

New

Ex vivo T cell CAR knock-in

RNA delivery



FL-1625T

DNA delivery

FL-1779T

New


Efficient IgG induction with saRNA in clinical trial (Ph 3 ongoing)

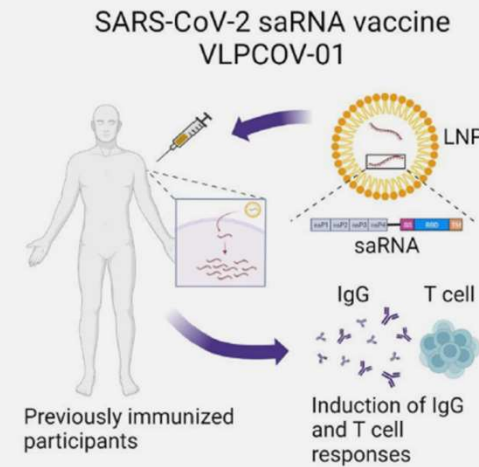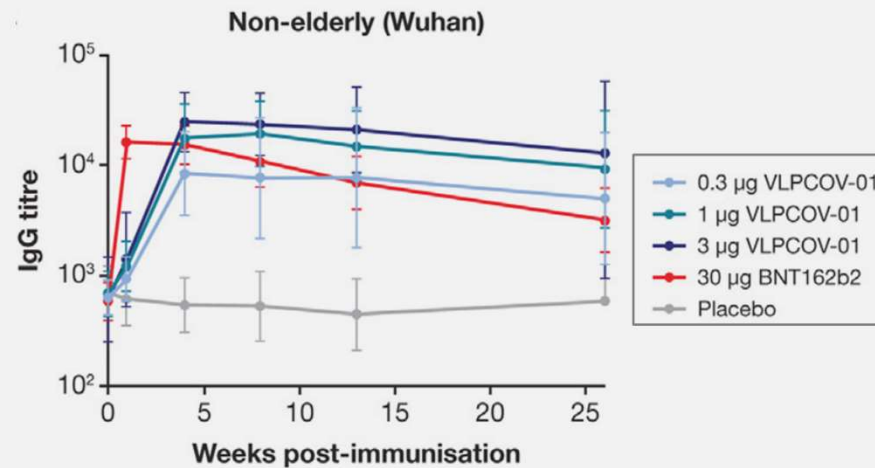

Cell Reports Medicine
2023, 4, 101134.

We completed process development and GMP Mfg. in both FL-0445 and its LNP

Ionizable lipid synthesis

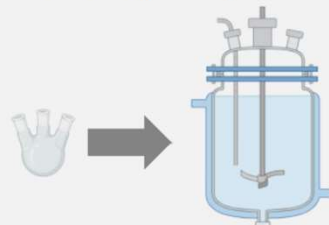

LNP production



FL-1245T | Tolerability test in rats

i.m. Vaccine	i.v. Liver	i.v. Liver- detarget.	Ex Vivo CAR-T
-----------------	-----------------------	-----------------------------	------------------

Poster 33

hEPO-mRNA



Rat (Sprague-Dawley, male, 7w, N=3)

3 mg/kg (mRNA), i.v. single dose

2023
FL-0207T

2024
FL-1207T

2025
FL-1245T

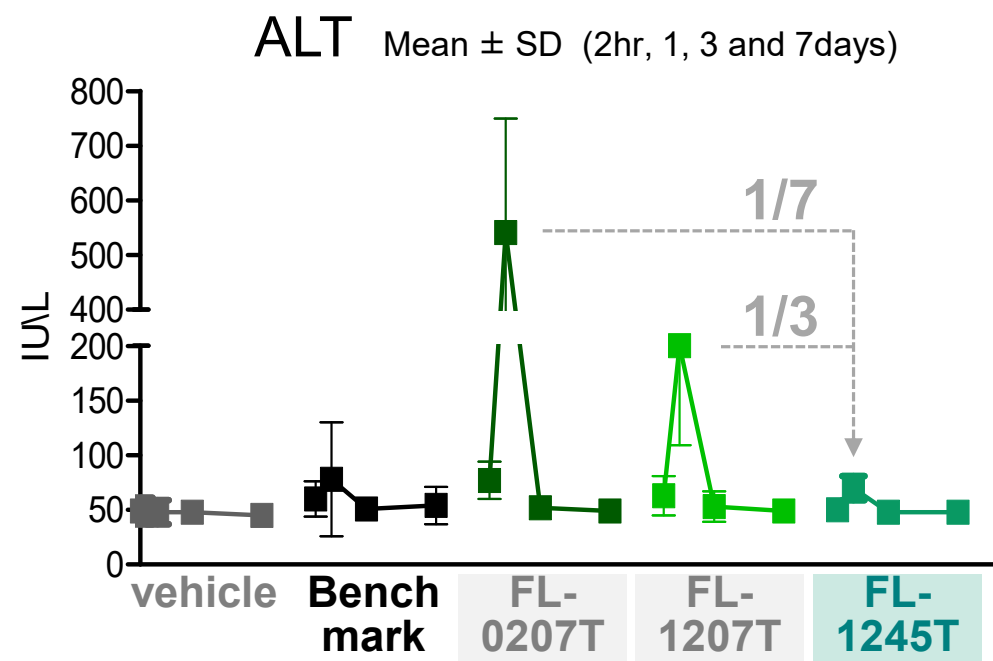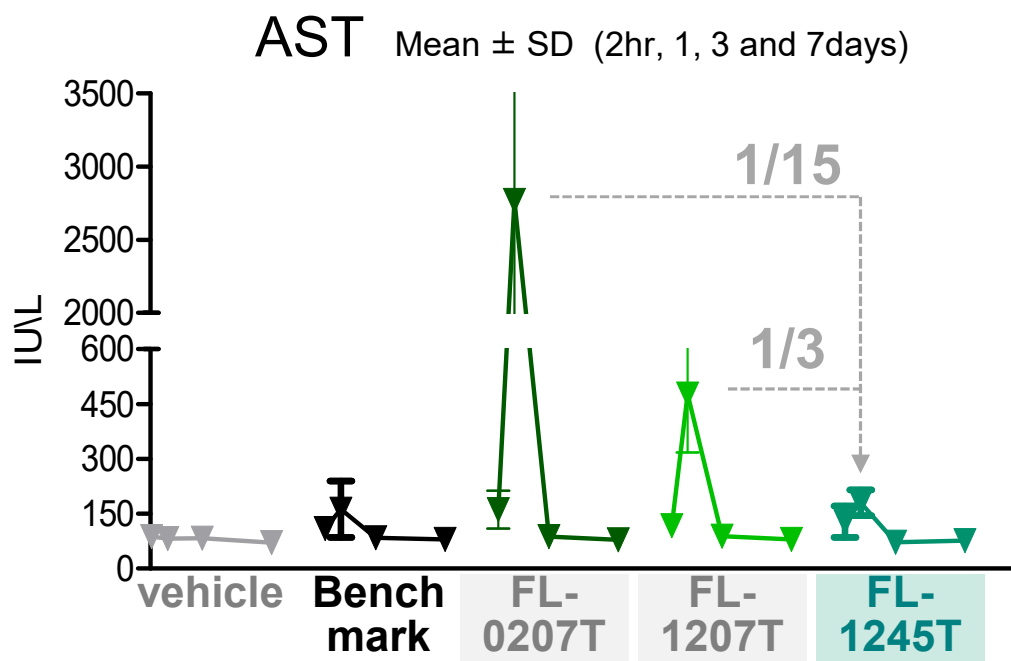

➤ **Our newest lipid FL-1245T demonstrated high tolerability**

FL-1245T | Tolerability test in rats

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.

Ex Vivo
CAR-T


Poster 33

hEPO-mRNA

Rat (Sprague-Dawley, male, 7w, N=3)

6 mg/kg (mRNA), i.v. single dose

Benchmark: 1/3 of the rats died within 1 day, **FL-1245T:** 1/3 of the rats died the next day.

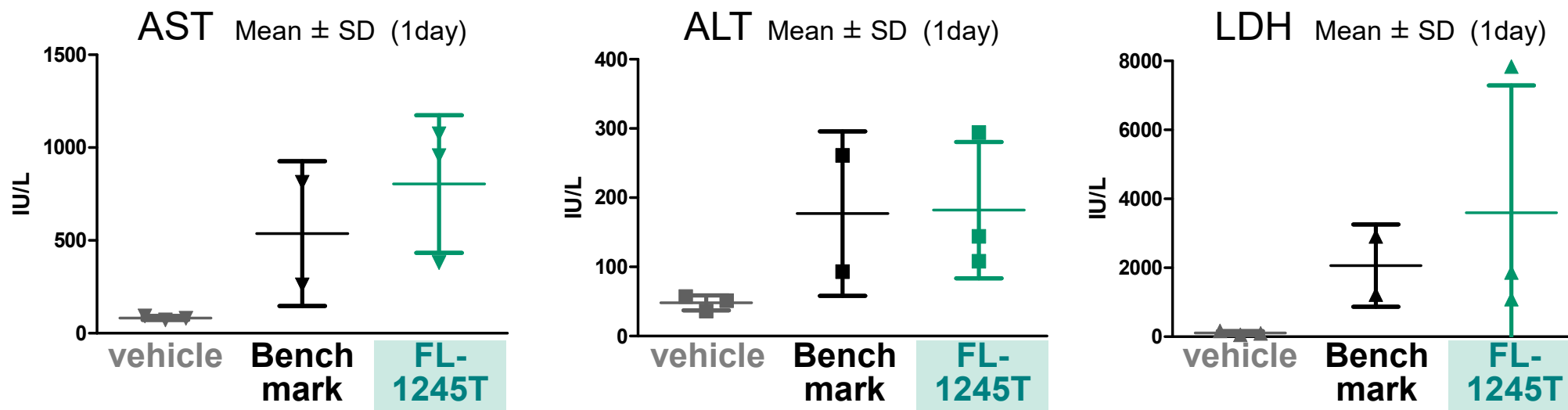

➤ **FL-1245T showed comparable tolerability to the benchmark lipid.**

FL-1245T | Protein expression in NHP



i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.


Ex Vivo
CAR-T

Poster 33

hEPO-mRNA

Cynomolgus monkey (male, 3y, N=3)

0.05 mg/kg (mRNA), 1hr infusion, i.v.

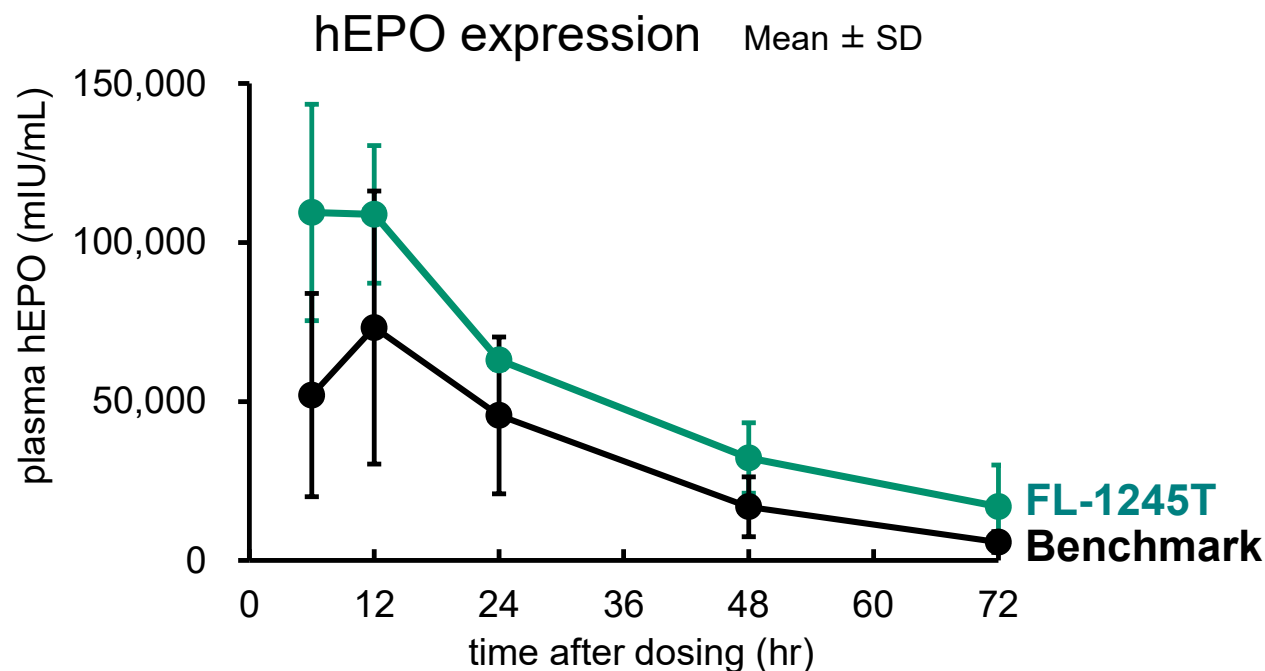

➤ **FL-1245T showed higher protein expression than benchmark lipid.**

We are planning higher dose studies


LNP potential in CAR-T cell manufacturing

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.

Ex Vivo
CAR-T

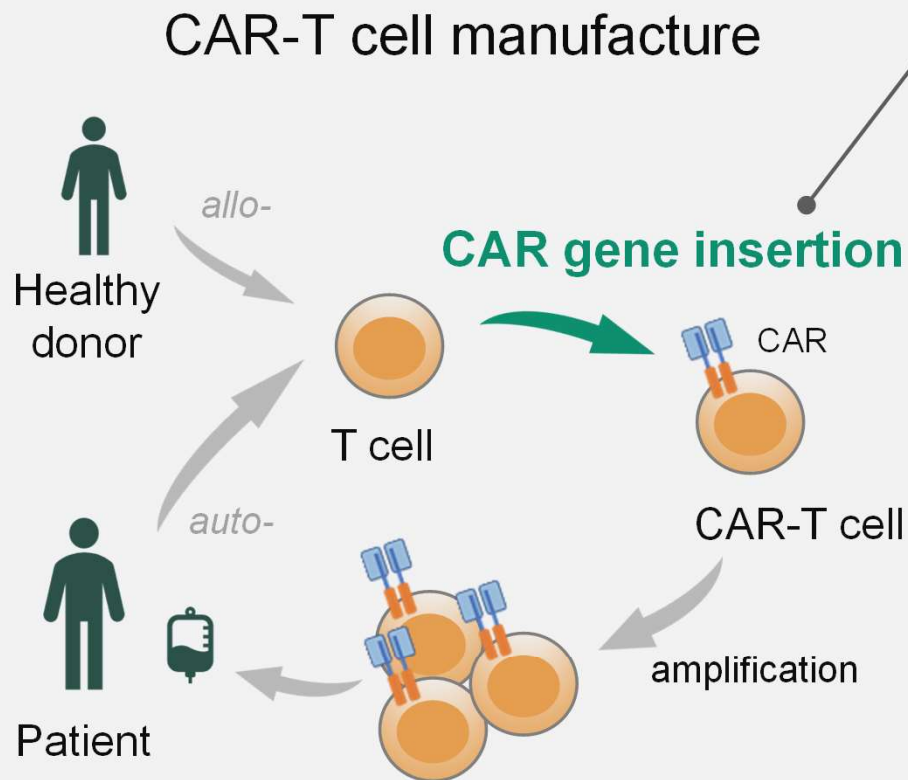


CAR gene insertion

Issues with current methods



- Viral production/QC high costs
- Safety concerns
- Cell damage by electroporation
- Low cell proliferation, low cell yield

Expectations for LNP methods

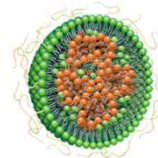

- No use of biological materials
- Reduce cytotoxicity, high yield

Pharmaceutics **2024**, 16(3), 346.

➤ **CAR gene insertion with LNP would overcome the current issues**



FL-1625T | T cell RNA delivery, knock-out

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.

Ex Vivo
CAR-T

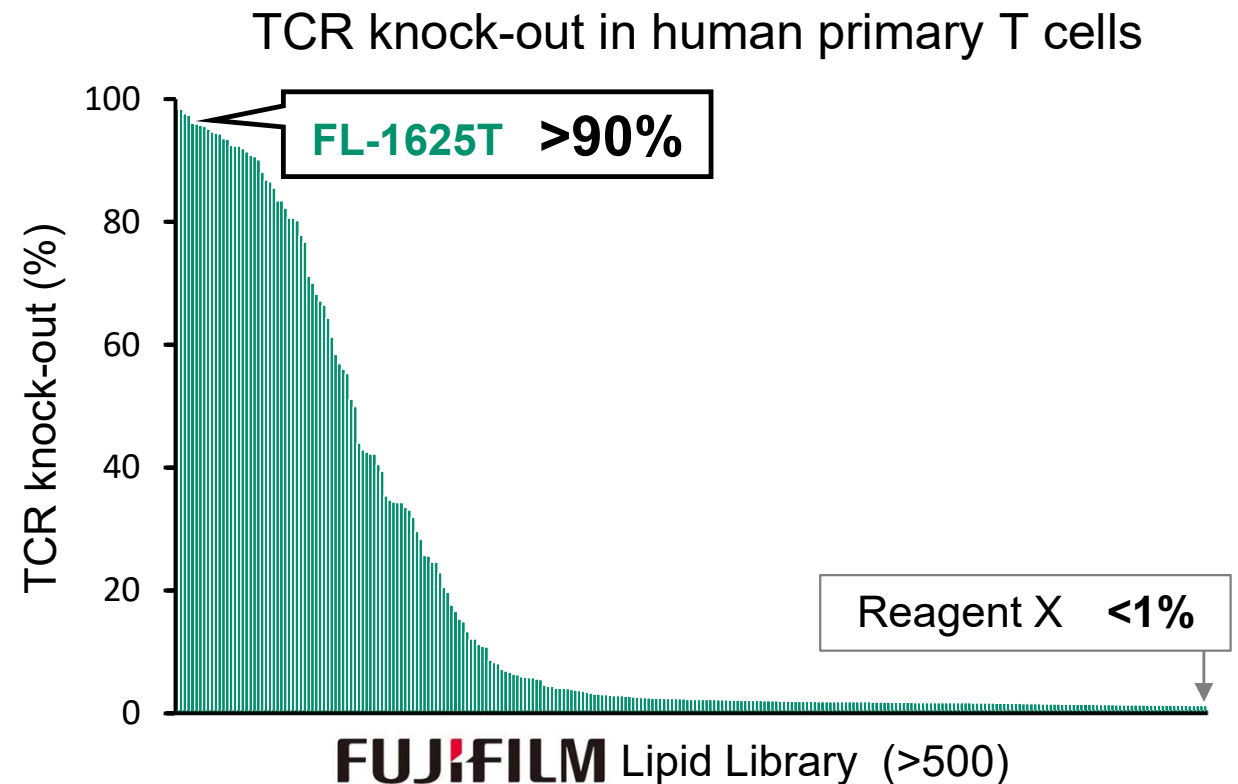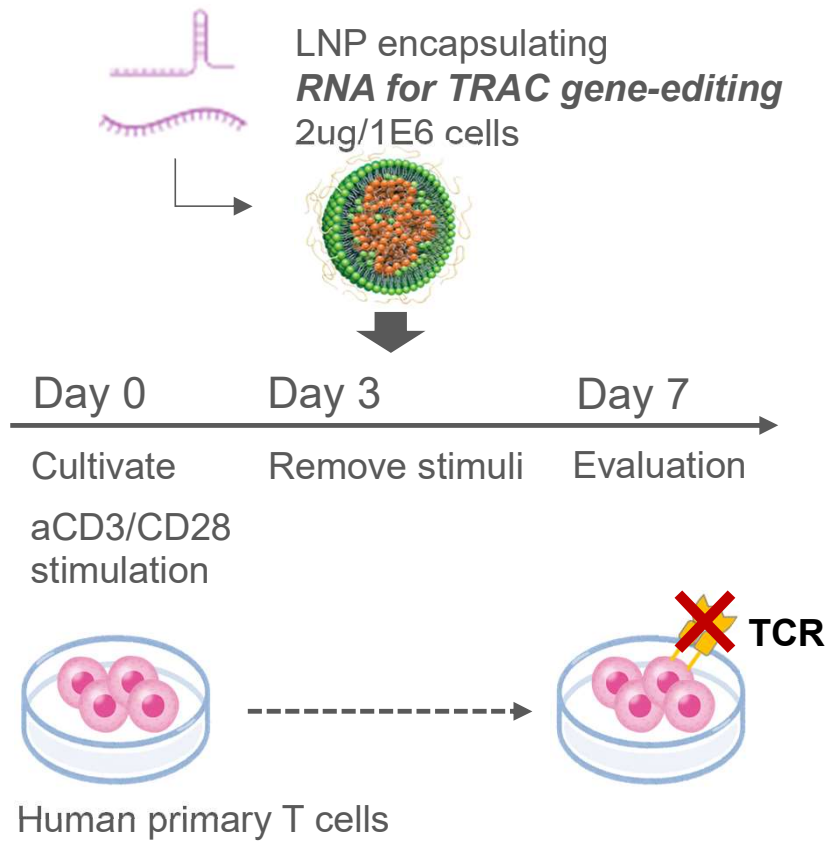

➤ **FL-1625T demonstrated high RNA delivery and >90% TCR knock-out.**



FL-1779T | T cell DNA delivery

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.

Ex Vivo
CAR-T

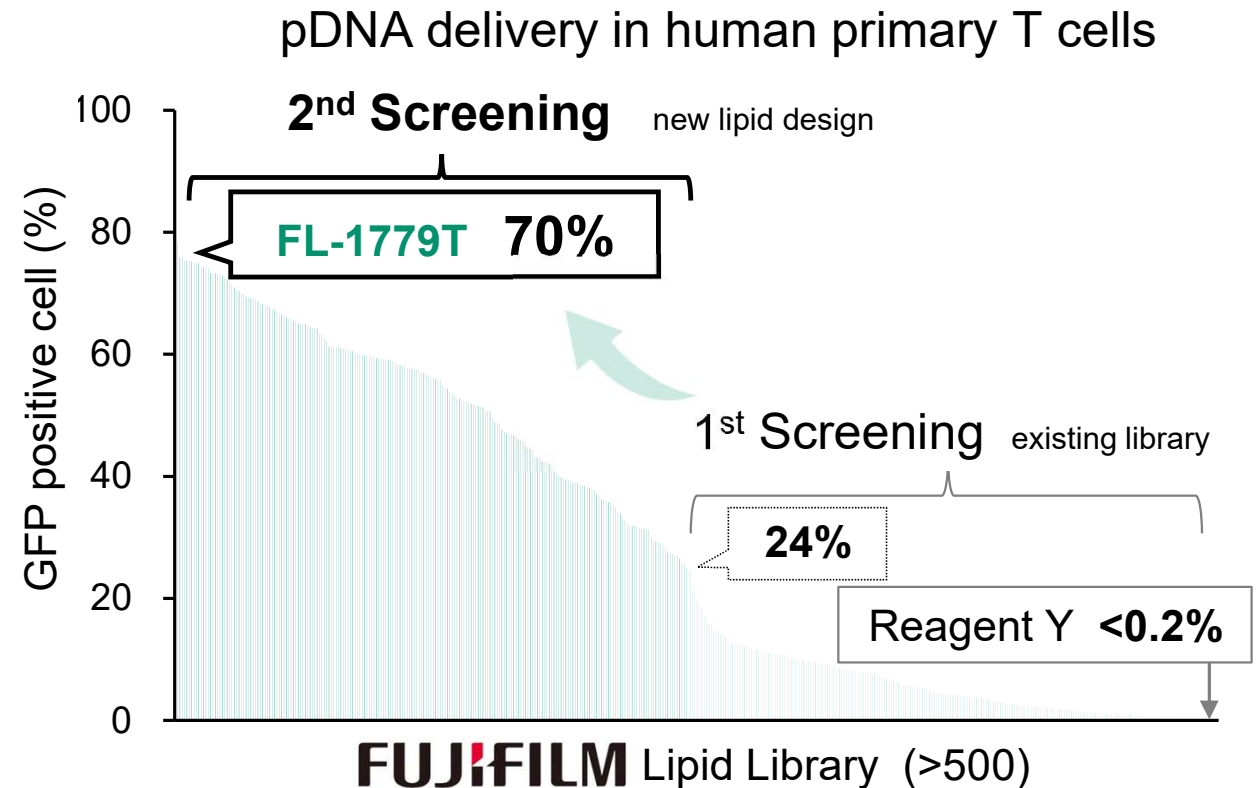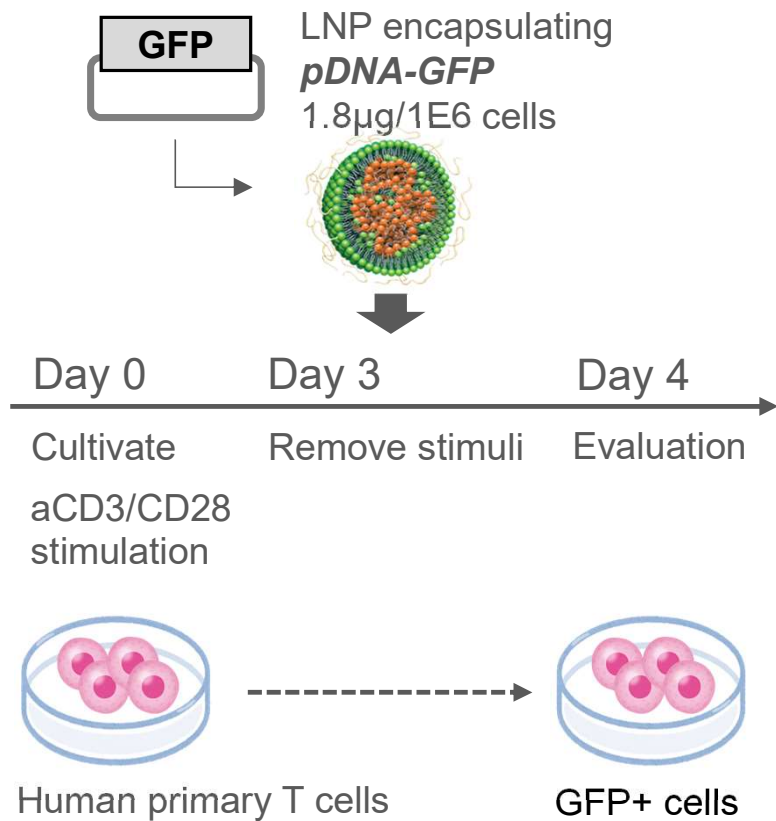

➤ We designed new lipid FL-1779T, which showed 70% pDNA delivery



FL-1625T×FL-1779T | T cell CAR knock-In

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.

Ex Vivo
CAR-T

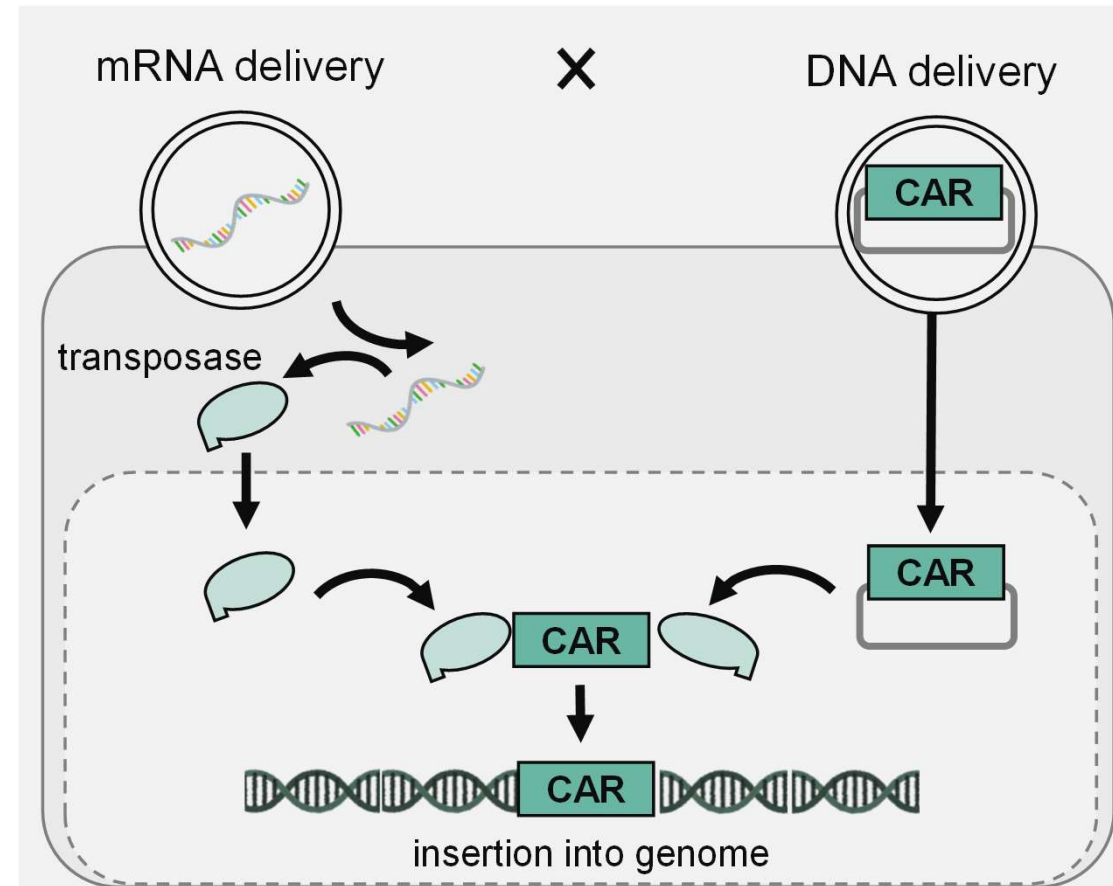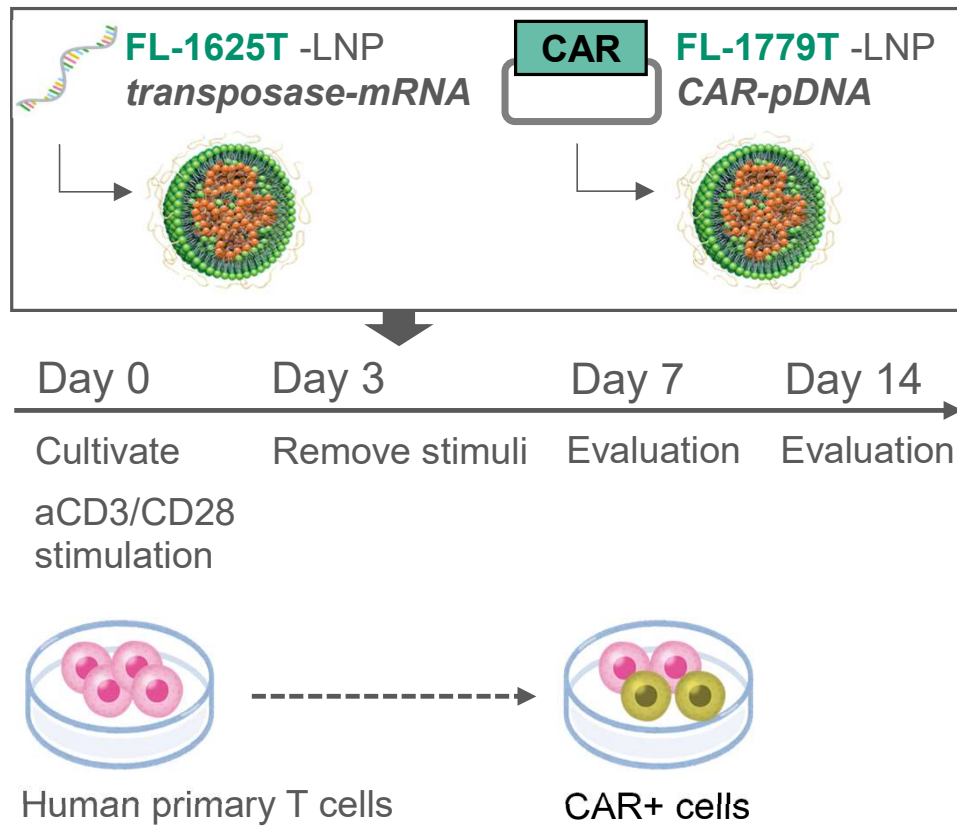

➤ By using two types of lipids, we tried CAR knock-in application.


FL-1625T×FL-1779T | T cell CAR knock-In

i.m.
Vaccine


i.v.
Liver

i.v.
Liver-
detarget.

Ex Vivo
CAR-T

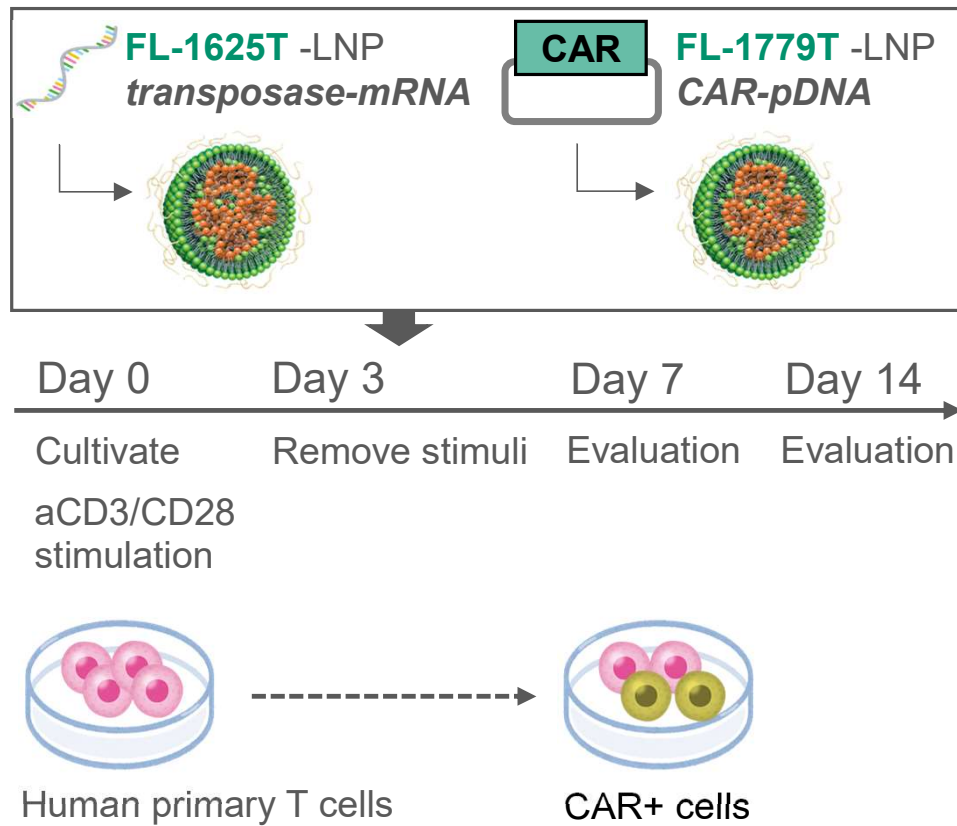

CAR-pDNA LNP/ transposase-mRNA LNP

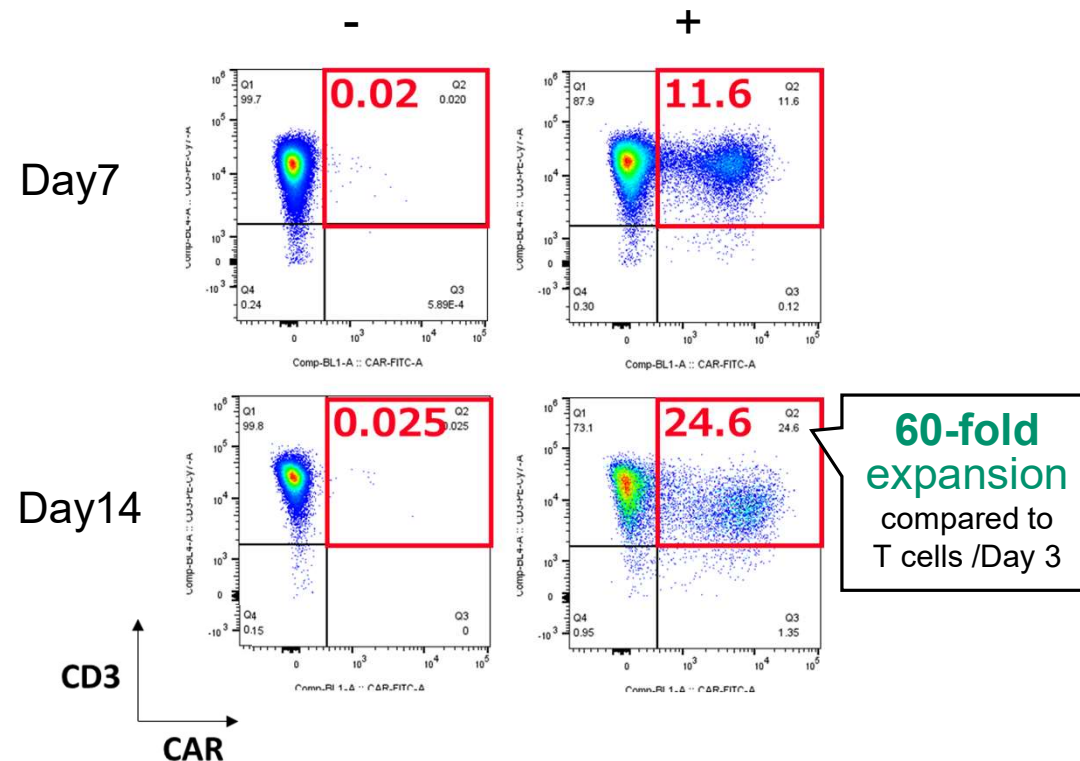

➤ CAR knock-in was achieved by using FL-1625T and FL-1779T.

FL-1010T | ApoE-independent uptake

i.m.
Vaccine

i.v.
Liver


i.v.
Liver-
detarget.

Ex Vivo
CAR-T


EGFP-mRNA

Ex vivo Activated T cell transfection

1.0μg/ 1E6 cells, ApoE (+/-)

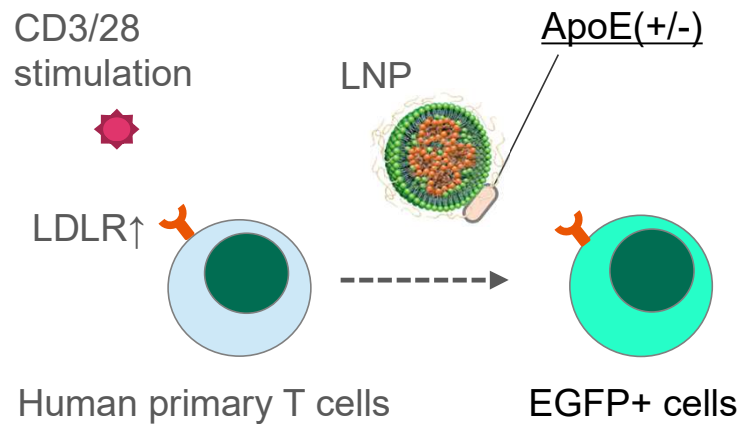

EGFP⁺ cell % in human primary T cells (1day)

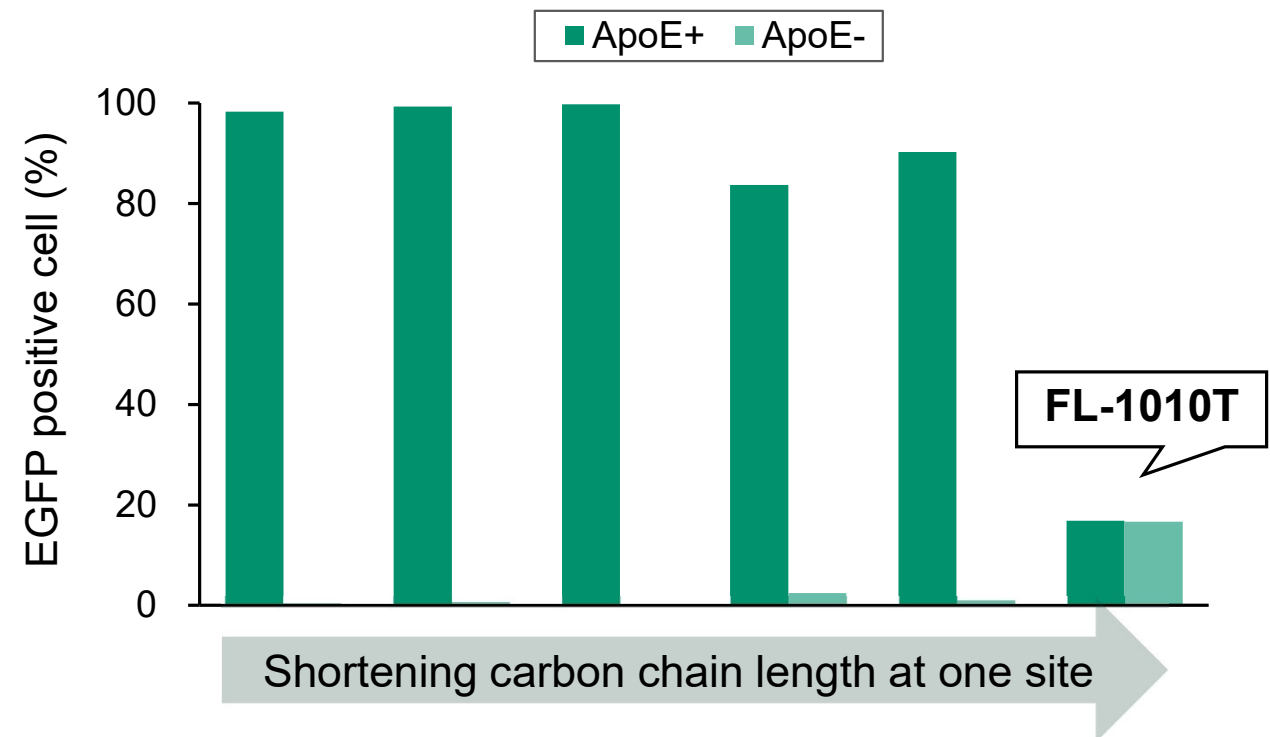

➤ **FL-1010T demonstrated ApoE independent uptake in ex vivo T cell study.**

Then, we conducted in vivo imaging experiments.

FL-1010T | In vivo expression distribution

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.


Ex Vivo
CAR-T

FLuc-mRNA

Mouse (ICR, female, 5wk, N=2)

0.2 mg/kg(mRNA), i.v.

Li: liver, Lu: lung, S: spleen K: kidney, H: heart
Circle Size indicates the approximate level of expression.
Expression Ratio: Based on Total Flux [p/s]

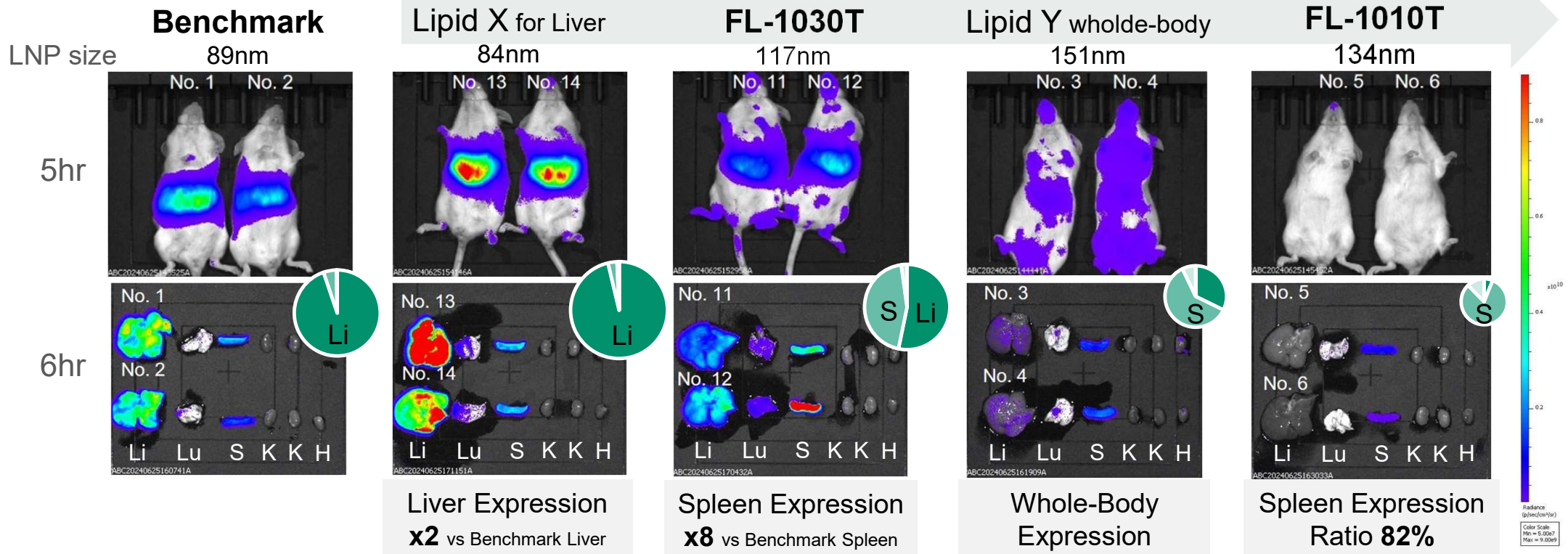

➤ Liver-detargeting was accomplished, but LNP size control was a bit delicate.

FL-1779T | ApoE-independent uptake

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.


Ex Vivo
CAR-T

EGFP-mRNA


Ex vivo Activated T cell transfection

1.0μg/ 1E6 cells, ApoE (+/-)

Poster 35

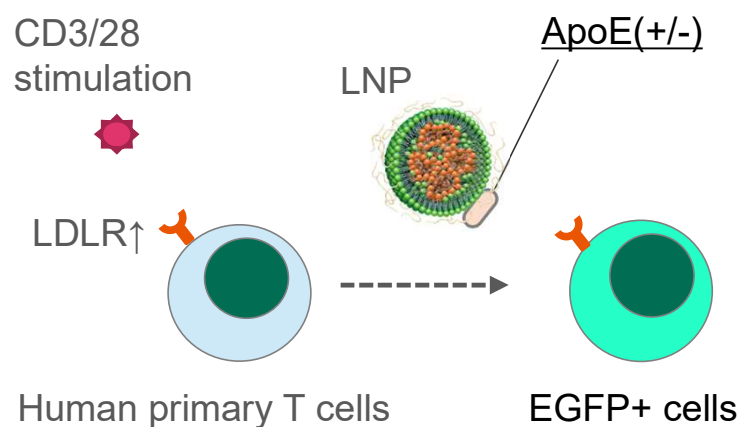

EGFP⁺ cell% in human primary T cells (1day)

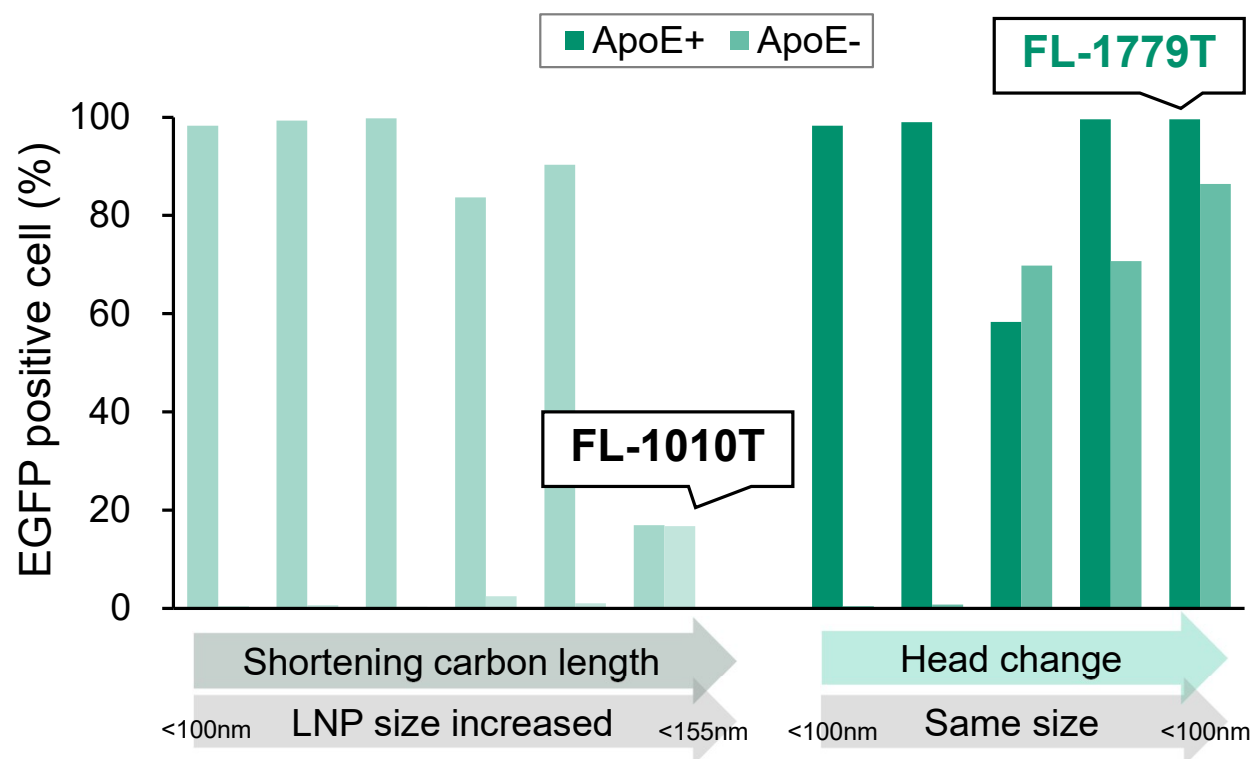

➤ **FL-1779T showed higher ApoE independent uptake and easier LNP size control.**

FL-1779T | Potential for extra-hepatic in vivo

i.m.
Vaccine

i.v.
Liver

i.v.
**Liver-
detarget.**

Ex Vivo
CAR-T



Poster 35

FLuc-mRNA

Mouse (ICR, female, 5wk, N=2)

0.2 mg/kg(mRNA), i.v. 6hr

Li: liver, Lu: lung, S: spleen K: kidney, H: heart

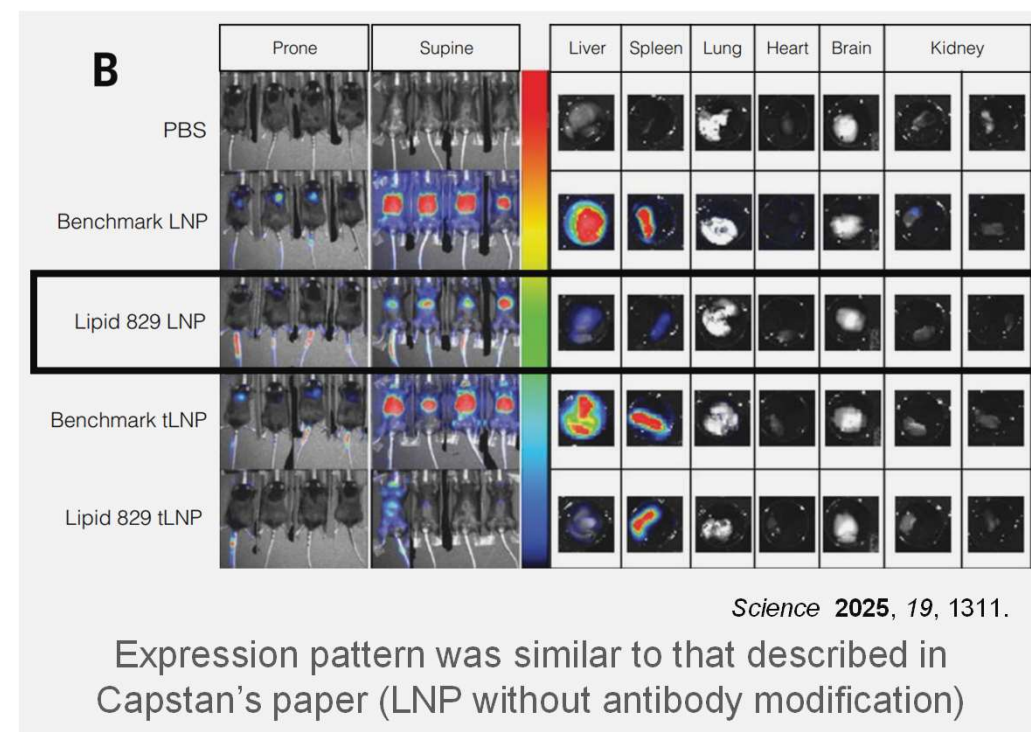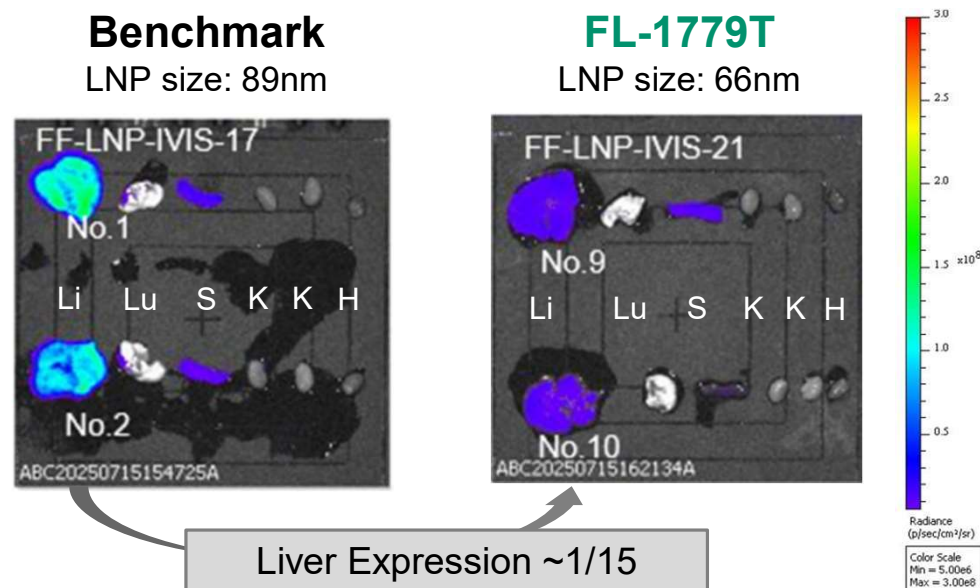

➤ **FL-1779T showed low expression in liver and spleen.**

Considering ex vivo data together, it may be suitable for in vivo T cell transfection (on-going)

FL-1779T×targeting-LNP | Ex vivo PBMC

i.m.
Vaccine

i.v.
Liver

i.v.
Liver-
detarget.

Ex Vivo
CAR-T


Poster 35

EGFP-mRNA


Specific IgG antibody-conjugated LNP

Ex vivo PBMCs transfection

ApoE (+)

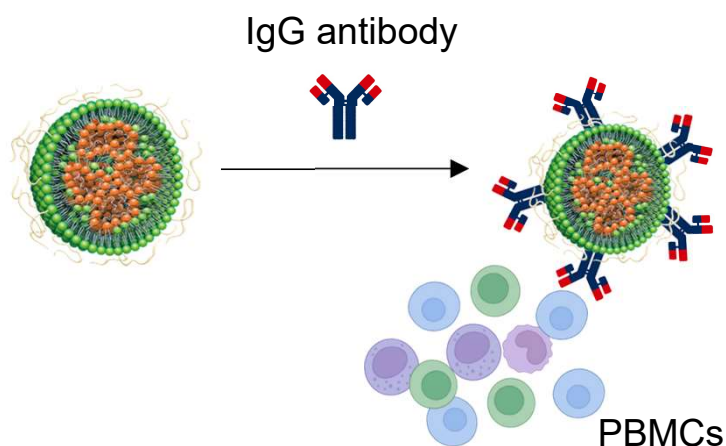

EGFP⁺ cell% in PBMC (1day)

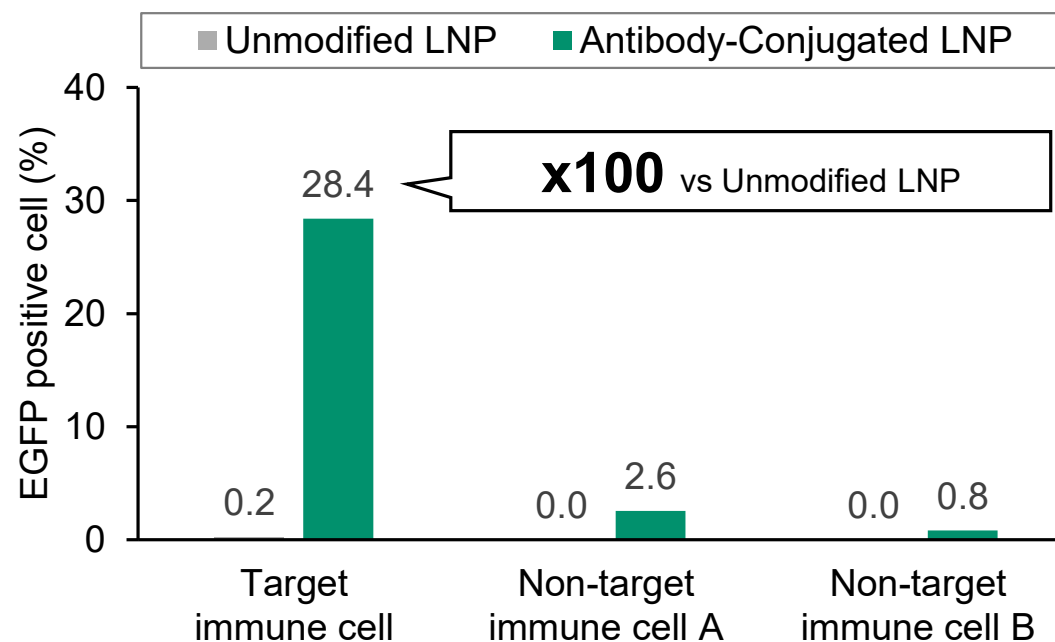

➤ FL-1779T-LNP conjugated with antibody showed immune cell selective expression.

Lipid Licensing

	Lipid	GMP Mfg.	IP	Remarks
i.m. vaccine	FL-0445	Established	Issued	Clinical trial (Ph 3)
	FL-2266	Established	Issued	GMP available
i.v. mRNA therapy	FL-1245T Poster 33		Pending	Liver mRNA delivery, high dose
	FL-1779T Poster 35		Pending	Liver de-targeting
	FL-1030T		Pending	Spleen mRNA delivery, high expression
Cell Therapy Gene Editing	FL-1625T		Pending	T cell RNA delivery (Empty LNP available)
	FL-1779T ---		Pending	T cell DNA delivery (Empty LNP available)

mRNA-LNP CDMO @Toyama/ Japan

Research	mRNA design and synthesis Poster 117 ~ mRNA-LNP formulation optimization
Development	process and analytical methods development
Manufacturing	mRNA-LNP manufacture in GMP



Acknowledgement

FUJIFILM Corporation

- Shigetomo Tsujihata
- Hirofumi Fukunaga
- Masaki Noro
- Yuta Murakami
- Nao Yamazaki
- Yuki Imaizumi
- Hirofumi Kimura
- Kentaro Numajiri
- Kohei Shimizu
- Kohei Yasuda
- Daisuke Nakagawa
- Sayako Umetani
- Misaki Masui
- Kazuhiko Nakata

and many colleagues

FUJIFILM Toyama Chemical Co., Ltd.

- Shinsuke Mizumoto
- Akira Inomata
- Hiromi Hayashi
- Keita Tanaka
- Tomomi Urabe
- Kaori Miyamoto
- Chie Kurosaki
- Takumi Koguchi
- Kazuki Himi
- Asuka Ejiri

and many colleagues

Fujifilm Group's Purpose

Giving our world more smiles

We bring diverse ideas, unique capabilities,
and extraordinary people together to change the world.

FUJIFILM
Value from Innovation

